Addendum

171/172 Diode Array Detectors User's Guide

Part Number LT350022

This document contains corrections and additions to information provided in the *171/172 Diode Array Detectors User's Guide*.

- The **Safety** and **Sécurité** sections replace the **Safety** chapter.

Safety

The instrument is intended to be used in a laboratory by trained technical personnel. For safe and correct use, both operating and service personnel should follow the instructions contained in the user's guide when installing, cleaning, and maintaining the instrument.

The following safety precautions must be observed during all phases of operation, service, and repair of the instrument. Failure to comply with these precautions or with specific warnings elsewhere in the user's guide violates safety standards of design, manufacture, and intended use of the instrument. Gilson assumes no liability for the customer's failure to comply with these requirements.

The instrument has been certified to safety standards required in Canada, Europe, and the United States. Refer to the rear panel label on the instrument or the Declaration of Conformity document for the current standards to which the instrument has been found compliant.

The following electronic and hazard symbols may appear on or inside the instrument.:

Symbol	Explanation	Symbol	Explanation
~	Alternating current Courant alternatif Wechselstrom		Caution Attention Vorsicht
	Protective conductor terminal Borne de terre de protection Schutzleiteranschluss		Caution, risk of electric shock Attention, risque de choc électrique Vorsicht, Elektroschockgefahr
	Electrical power ON Sous tension Netzschalter ein		Fuse Fusible Sicherung
o	Electrical power OFF Hors tension Netzschalter aus		Caution, hot surface Attention, surface chaude Vorsicht, heiße Oberfläche
	Direct current Courant continu Gleichstrom		Caution, ultraviolet light Attention, rayonnement ultraviolet Vorsicht, Ultraviolettes Licht

Voltage

Access to the rear panel is necessary. The instrument must be detached from all voltage sources before service, repair, or exchange of parts. For normal operation, the instrument is to be grounded through the AC line cord and power supply provided. Failure to do so can result in a potential shock hazard that could result in serious personal injury.

Use only fuses with the rated current and of the specified type as listed on the rear panel label on the instrument. The instrument must only be operated with the voltage specified on the rear panel label of the instrument and with the grounded AC line cord and power supply provided.

Heat

The deuterium lamp can reach temperatures of 176°F (80°C) in normal operation. Allow 30 minutes for the lamp to cool before servicing/replacing.

Solvents

Observe safe laboratory practices when handling solvents. If working with hazardous solvents or flammable liquids, ensure that there is proper ventilation and that adequate protection such as safety glasses, gloves, and protective clothing are used.

If dangerous liquids are used, precautions should be taken to limit potential hazards from leaks and/or spillage through the use of a non-flammable tray or use of a fume hood, etc.

If there is the potential of explosive gases being developed, a fume hood or other means should be used to safely manage that risk.

Refer to the Material Safety Data Sheets for the solvents before use.


Replacement Parts

Be sure to use only replacement parts mentioned in the user's guide. Do not repair the instrument or change parts not listed in the user's guide. If it is necessary to replace parts not listed, please contact your local Gilson representative.

Ultraviolet (UV) Radiation

WARNING Never look directly into the light from the deuterium lamp. Always use protective glasses with a UV filter, if you have to open the lamp housing while the deuterium lamp is on.

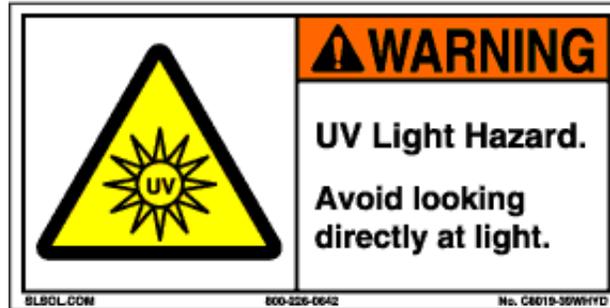Warning: The deuterium lamp emits UV radiation that is harmful to the eyes. Always turn off the detector power before removing the lamp cover.

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Fiber Optics

WARNING Do not look directly at the light emitted from the fiber optics.

Please note the following when handling the fiber optics:

- Observe a minimum bend radius for optical fibers. For a fiber optic with a diameter of 0.6 mm, the smallest allowable bending radius is approximately 50 mm for internal fibers and approximately 120 mm for external fibers.
- Avoid contact with fingers or contamination of the optical fiber tip. Even very thin layers of contaminants will greatly impair measurements, particularly in the UV range.
- If the front surface of the optical fiber tip becomes dirty, it should be cleaned with a lint-free cloth.

When not in use, the ends of the optical fibers should always be protected with the end caps delivered with the fibers.

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