Quick Guide

EN | FR

Parts Checklist

	COMPONENT	MATERIAL
1	Aspiration hand operator	Delrin® acetal resin and silicon
not pictured	Silicon adapter for Pasteur pipette (already mounted under the bayonet cap)	Silicon
2	Safety stand for the hand operator with storing gaps for the adapters	Delrin acetal resin
3	Single-channel plastic tip adapter with ejector	Delrin acetal resin
4	8-channel plastic tip adapter with ejector	Delrin acetal resin
5	8-channel stainless steel needle adapter	Delrin acetal resin and stainless steel
6	Four single-channel stainless steel needle adapters (various sizes): <ul style="list-style-type: none"> • 4 cm length, 1.5 mm diameter • 4 cm length, 3 mm diameter • 12 cm length, 1.5 mm diameter • 12 cm length, 3 mm diameter 	Delrin acetal resin and stainless steel
not pictured	Silicon adapter for Pasteur pipette	Silicon
not pictured	Silicon grease	Silicon

Image 1

Safe Aspiration Kit components
Composants du Safe Aspiration Kit

Image 2

Installation steps
Etapes d'installation

Installation

- A** Insert the back of the aspiration hand operator with barbed fitting into the silicon tubing of the aspiration system. (Refer to Image 2)
- B** Insert a disposable aspiration pipette, the single- or 8-channel stainless steel adapter directly into the port on the front of the aspiration hand operator. (Refer to Image 2)
For the single- or 8-channel plastic tip adapter with ejector, remove the bayonet cap and silicon adapter. Then screw on the plastic tip adapter with ejector. (Refer to Image 2)
- C** Vacuum pressure can be controlled by depressing dark blue button on top of the aspiration hand operator. (Refer to Image 2)

NOTICE Residual vacuum can be controlled by turning dark blue plastic screw at bottom of the aspiration hand operator. Turning screw one full turn clockwise shuts the valve and cuts off the vacuum. Turning screw counter-clockwise adjusts the constant suction level.

NOTE The safe aspiration hand operator is factory pre-set to allow a low level of constant suction.

Cleaning and Special Care

- The safe aspiration hand operator and accessories are fully autoclavable.
- Clean by immersing all pieces in a solution of water and mild laboratory detergent. Do not use abrasives, as this may affect operation of the control valve. Rinse with distilled water.
- Autoclave the unit prior to disassembly for cleaning after working with biohazardous materials. Refer to Image 3 for disassembly instructions.
- Immediately flush the aspiration hand operator with ethanol after working with organic solvents.
- To avoid build-up of deposits, immediately flush the aspiration hand operator with distilled water after working with salt-rich solutions.
- Apply a thin coat of silicon grease as needed to the control valve. (Refer to Figure 3)

Recommended Tips

For optimum performance of the single- and 8-channel plastic tip adapters with ejector, we recommend using PIPETMAN® DIAMOND Tips DL10, D200, and D300.

Warranty

Gilson warrants this Safe Aspiration Kit against defects in material under normal use and service for 1 year from the date of purchase. This warranty is valid only if the components of the kit are used in the manner described in this guide and for the purpose for which it is designed. Gilson is not responsible for consequential damages resulting from the misuse or bad cleaning or decontamination of this instrument. Forced opening of the instrument invalidates any warranty claim. Enclose with the returned instrument a description of the trouble that has occurred and specify which reagents were used.

NOTICE Please clean the instrument carefully before returning it. Never send in instruments filled with reagents. Returned instruments can only be checked and repaired if they have been carefully cleaned and decontaminated by the customer. Note that freight costs related to the return of the instrument will be the responsibility of the customer.

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Specifications subject to change without notification — errors excepted. Reproduction, adaptation, or translation without prior written permission is prohibited, except as allowed under copyright laws.
LT801604/A | ©2018 Gilson, Inc. All rights reserved.