Quick Guide

EN | FR

*Discover the new experience
of connected pipetting on
www.gilsonconnect.com*

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Parts Checklist

- SAMSUNG Galaxy Tab S3 with PipettePilot application
- Power adapter, 5V
- Tempered glass screen protector
- Tablet holder
- Microplate holder
- Microplate adapter
- Strip adapter
- Environment sensor
- TRACKMAN® Connected Quick Guide

Getting Started

NOTICE

Charge the SAMSUNG Galaxy Tab S3 before first use.

- 1 Insert the smaller part of the tablet holder, into one of the five slots of the larger part according to the inclination you need. According to your personal comfort, you can choose one of the two tablet holder positions below.

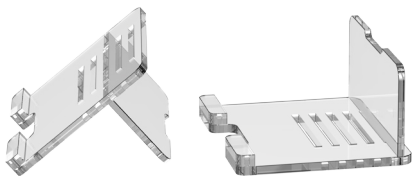
Figure 1

Both options of tablet holder assembly

- 2 Place the SAMSUNG Galaxy Tab S3 on the tablet holder and turn it on.

NOTICE

It is strongly recommended to install the screen protector provided before use

NOTE

The tablet user manual is available directly on the tablet.

User manual
User manual

- 3 Launch PipettePilot application.
- 4 Turn the environment sensor on by pressing gently on the purple side. The LED will blink three times to indicate that it is on.

When prompted by PipettePilot, connect the environment sensor to the tablet using the Bluetooth® connection. Once turned on and connected, the environment sensor will automatically begin monitoring temperature, humidity, and pressure.

To turn off the environment sensor, press and hold the purple side down for five seconds. The LED will blink rapidly to indicate that it is turned off.

NOTE

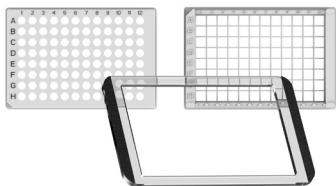
The environment sensor user manual is available at www.bluemaestro.com/tempo-disc-temperature-humidity-resources/

NOTICE

Pressing the button with excessive force may damage the device.

**Figure 2**

Environment sensor

**Figure 3**

Microplate holder (middle), microplate adapter (left), and strip adapter (right)

- 5 When prompted by PipettePilot, remove the two white strips from the microplate holder and then place the microplate holder on the corresponding frame on the tablet. If needed, use the microplate adapter or the strip adapter.
- 6 Follow PipettePilot instructions.

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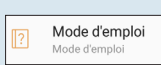
Other References

This Quick Start Guide provides you with basic information to get started with TRACKMAN Connected.

Make sure you have read the TRACKMAN Connected User's Guide, available at: www.gilsonconnect.com and other user manuals provided with the environment sensor (available at www.bluemaestro.com/tempo-disc-temperature-humidity-resources/), screen protector, and tablet (inside the box).

Warranty

Gilson warrants the devices against defects in material under normal use and service for a period of 12 months from the date of purchase.

This warranty shall not apply to devices which are subject to abnormal use and/or improper or inadequate maintenance (contrary to the recommendations given in the user's guide), including, but not limited to device which have been subjected to physical damage, improper handling, or spillage or exposure to any corrosive environment. This warranty shall also be void in the event devices are altered or modified by any party other than its designates. Gilson's sole liability under this warranty shall be limited to, at Gilson's sole option, repair or replacement of any defective components of the devices or refund of the purchase price paid for such device.

THE FOREGOING WARRANTY IS EXCLUSIVE AND GILSON HEREBY DISCLAIMS ALL OTHER WARRANTIES, WHETHER EXPRESS OR IMPLIED, INCLUDING ANY WARRANTIES OF MERCHANTABILITY AND ANY WARRANTIES OF FITNESS FOR A PARTICULAR PURPOSE, UNDER NO CIRCUMSTANCES SHALL GILSON BE LIABLE FOR ANY CONSEQUENTIAL, PUNITIVE, INDIRECT OR INCIDENTAL DAMAGES ARISING OUT OF ANY BREACH OF ANY EXPRESS OR IMPLIED WARRANTY.

Warranty for Non-Gilson Equipment

Contact your Gilson representative. Gilson will coordinate with the manufacturer for handling warranty-related service. The warranties are those provided by the manufacturer.

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LT801593/A | ©2018 Gilson. All rights reserved.